Look for these other great products from Learning Resources®:

LER 1963 Easy Grip Tweezers (Set of 12)

LER 2784 Primary Science Set

www.LearningResources.com

Visit our website to write a product review
or to find a store near you.

CE © Learning Resources, Inc., Vernon Hills, IL (U.S.A.)
Learning Resources Ltd., King's Lynn, Norfolk (U.K.)
Please retain our address for future reference.
Made in China. LRM2933-GUD

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3 Way Magnification Science Station

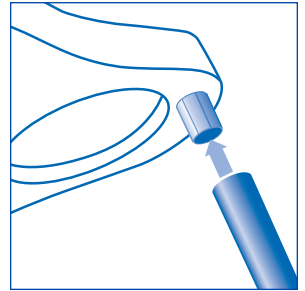
Contents

- Lens Frame with (3) Acrylic Lenses
- 3 Removable Legs

Assembly

Gently push the legs onto the posts on the bottom of the frame. You may need to twist the legs slightly. Rubber caps on the legs help the frame remain steady and limit skidding.

To remove the legs, gently twist and pull them off the frame.

Suggested Use

The 3-Way Magnification Science Station is a great tool for hands-on exploration. The varying magnifications allow for easy inspection of objects of all shapes, sizes, and surfaces.

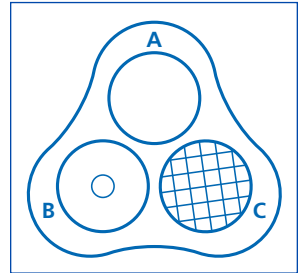
Simply place an object underneath the lens, on a table or flat surface, and view from overhead. Depending on its size and shape, the object may need to be lifted up until it comes into focus.

A. 4X Magnifying Lens

This lens is best suited for exploring general objects.

B. 8X Dual-Magnifying Lens with 20X Super Magnifier

Use this lens to compare and contrast magnifications and to focus in closely on objects with fine detail. Note: You may need to adjust the distance between the object and the inner 20X lens because of its strong magnification. Try closing one eye to help focus.

C. Distortion Lens w/1X Magnification

This lens mimics the compound eye of many insects. Objects will be seen as multiple images. Try twisting and turning objects while underneath the lens to increase the effect.

General Care

To clean lenses, always use a clean, soft cloth. Avoid using soap or solvents that can build up and cloud the lenses.

The lenses in this product are made of acrylic. Please take precautions to avoid dropping or scratching the lenses.

Caution: Never use magnification lenses to look at the sun or other light sources that can damage vision.

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